MC12

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Flexicon
Liquid Filling

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
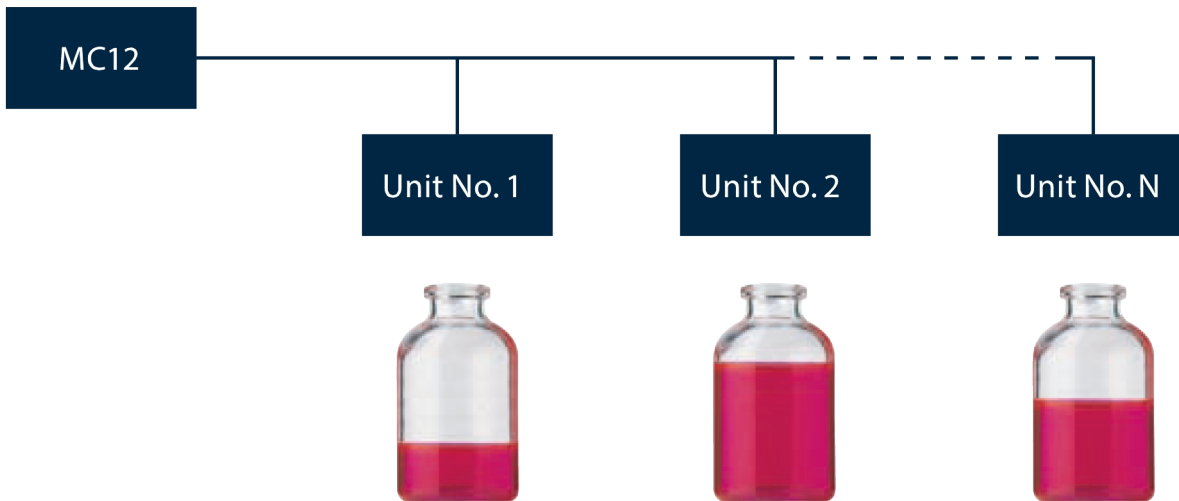
INDIVIDUAL MODE

MC12 can control up to 16 fillers regardless of the type. MC12 controls each filler using individual parameters such as fill volume, tube

diameter, acceleration, use of a sensor for »no bottle – no fill«, activation by an external signal and more. Each filler is calibrated individually.

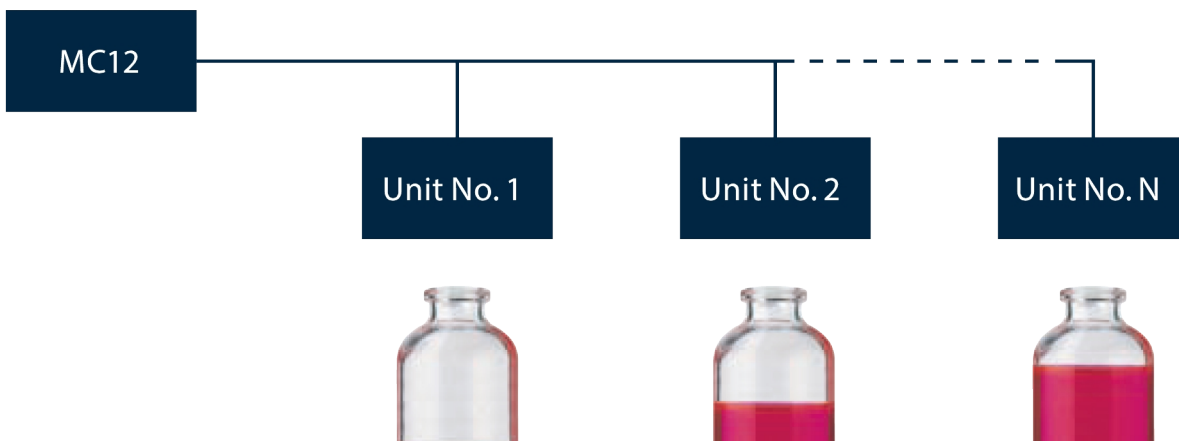
SERIAL MODE

In serial mode, the total fill is made up from the accumulated volumes delivered by each filling station. The fillers can operate with individual tube diameters, thus giving different fill volumes from each filling needle.

Calibration is carried out on all filling stations simultaneously. A sensor for »no bottle – no fill« may be connected. This set-up is used for medium to large production in fully automatic filling systems, primarily using a star wheel.

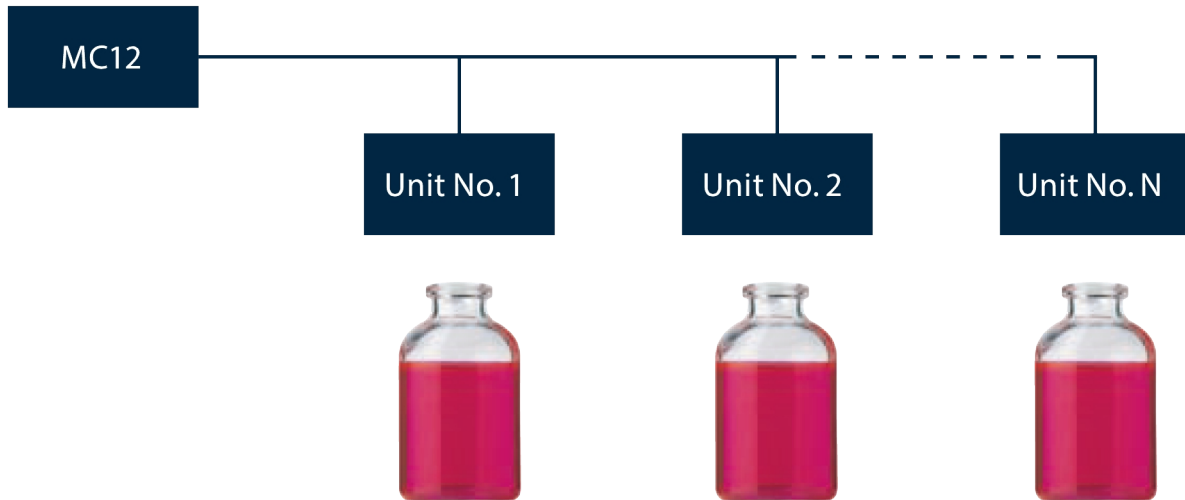PARALLEL MODE

In parallel mode, all connected fillers are synchronized with the same set of parameters. The calibration is carried out individually.

A sensor for »no bottle – no fill« may be connected. This set-up is used for larger production volumes, primarily filled on in-line filling systems.

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